


# Extraction of Synthetic Cannabinoids (SPICE) from Whole Blood using ISOLUTE® SLE+ prior to GC-MS Analysis

This application note describes the extraction of a range of synthetic cannabinoids and metabolites from whole blood prior to GC-MS analysis. An effective and efficient ISOLUTE® SLE+ protocol has been developed that is optimized for extraction of 800 µL of pre-treated matrix. The simple sample preparation procedure delivers clean extracts and analyte recoveries greater than 85% with RSDs lower than 10% for all analytes.

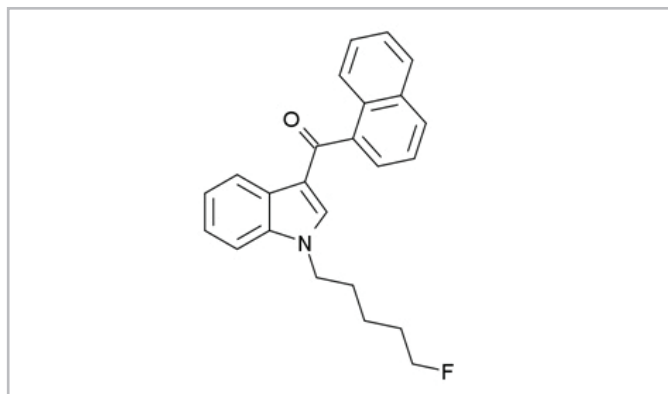

**Figure 1.** An example of a synthetic cannabinoid, AM-2201

## Introduction

ISOLUTE SLE+ Supported Liquid Extraction plates and columns offer an efficient alternative to traditional liquid-liquid extraction (LLE) for bioanalytical sample preparation, providing high analyte recoveries, no emulsion formation, and significantly reduced sample preparation.

## Analytes

UR-144, JWH-073, JWH-018, 5-hydroxypentyl-JWH250, 3-hydroxybutyl-JWH073, AM-2201, 4-hydroxypentyl-JWH-018, 5-hydroxypentyl-JWH-018, JWH-200

## Sample Preparation Procedure

**Sample Pre-treatment** Dilute whole blood (400 µL) with water (400 µL) and vortex for 10 seconds.

### ISOLUTE SLE+ 1 mL Sample Volume columns, part number 820-0140-C

**Sample Loading:** Load the pre-treated whole blood (800 µL total volume) onto the column and apply a pulse of vacuum or positive pressure to initiate flow. Allow the sample to adsorb for 5 minutes.

**Analyte Extraction:** Apply hexane/ethyl acetate (95/5, v/v, 2.5 mL) and allow to flow under gravity for 5 minutes. Apply a further aliquot of hexane/ethyl acetate (95/5, v/v, 2.5 mL) and allow to flow for another 5 minutes under gravity. Apply vacuum or positive pressure to pull through any remaining extraction solvent.

**Post Elution and Derivatisation:** Evaporate to dryness in a stream of air or nitrogen using a SPE Dry (40 °C, 20 to 40 L/min) or TurboVap (1.5 bar at 40 °C for 40 mins). Reconstitute with ethyl acetate (250 µL) and vortex for 20 seconds. Transfer to a high recovery glass vial and evaporate to dryness.

Add ethyl acetate (25 µL) and BSTFA:TMCS 99:01 (25 µL) and cap with a non-split cap. Vortex for 20 seconds and heat in a heating block set to 70 °C, for 30 minutes. Remove vial from the block and allow to cool.

## GC Conditions

<b>Instrument:</b>	Agilent 7890A with QuickSwap
<b>Column:</b>	SGE capillary column; BPX5, 30 m x 0.25 mm ID x 0.25 µm
<b>Carrier:</b>	Helium 1.2 mL/min (constant flow)
<b>Inlet:</b>	250 °C, Splitless, purge flow: 50 mL/min at 1.5 min, septum purge flow: 3 mL/min
<b>Injection:</b>	2 µL
<b>Wash Solvent:</b>	Ethyl acetate
<b>Oven:</b>	Initial Temperature 100 °C Ramp 50 °C/min to 275 °C, hold for 4 minutes Ramp 120 °C/min to 300 °C, hold for 8 minutes Ramp 120 °C/min to 315 °C, hold for 1.5 minutes Ramp 120 °C/min to 330 °C, hold for 1.5 minutes
<b>Post Run:</b>	Backflush for 2.4 minutes (3 void volumes)
<b>Transfer Line:</b>	280 °C

## MS Conditions

<b>Instrument:</b>	Agilent 5975C
<b>Source:</b>	230 °C
<b>Quadrupole:</b>	150 °C
<b>MSD mode:</b>	SIM


## SIM Parameters

**Table 1.** Ions acquired in the Selected Ion Monitoring (SIM) mode


SIM Group	Analyte	Target (Quant) Ion	1 <sup>st</sup> Qual Ion	2 <sup>nd</sup> Qual Ion	3 <sup>rd</sup> Qual Ion
1	UR144	214	311	N/A	N/A
2	RCS-4	320	263	N/A	N/A
3	JWH073	327	200	310	N/A
4	JWH018	341	214	324	N/A
5	5-hydroxypentyl-JWH250	302	228	N/A	N/A
6	3-hydroxybutyl-JWH073	285	270	415	N/A
7	AM2201	359	284	342	N/A
8	4-hydroxypentyl-JWH018	284	270	296	429
9	5-hydroxypentyl-JWH018	270	284	414	N/A
10	JWH200	100	384	N/A	N/A

## Results


This optimized ISOLUTE® SLE+ protocol demonstrated analyte recoveries ranging from 85-99% from whole blood, as shown in **Figure 2**. RSDs were below 10% for all analytes.

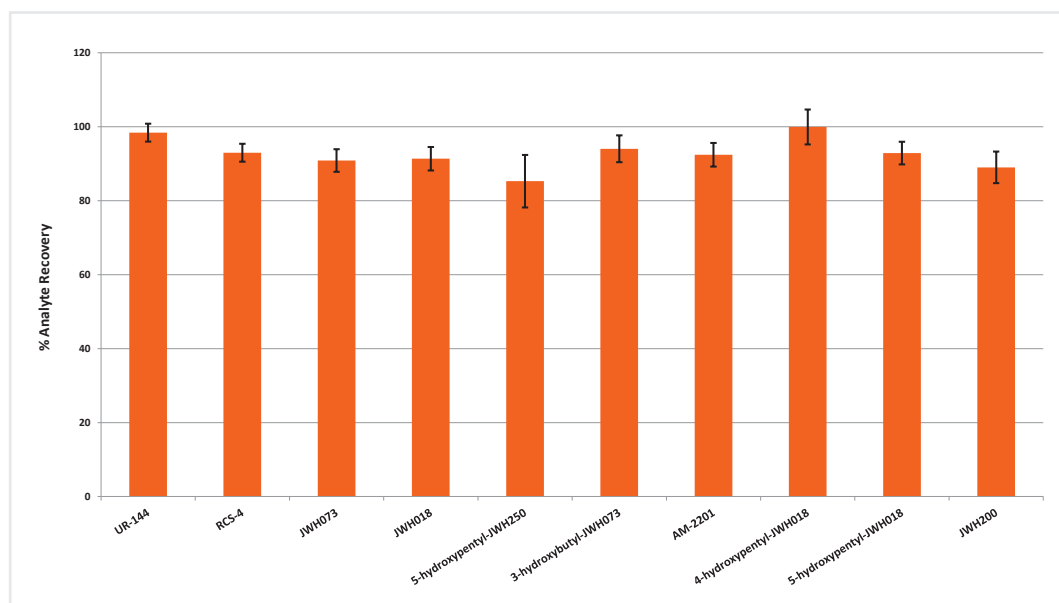

**Figure 2.** Typical analyte % recoveries for extracted synthetic cannabinoids and metabolites (n=7) using the ISOLUTE® SLE+ protocol.

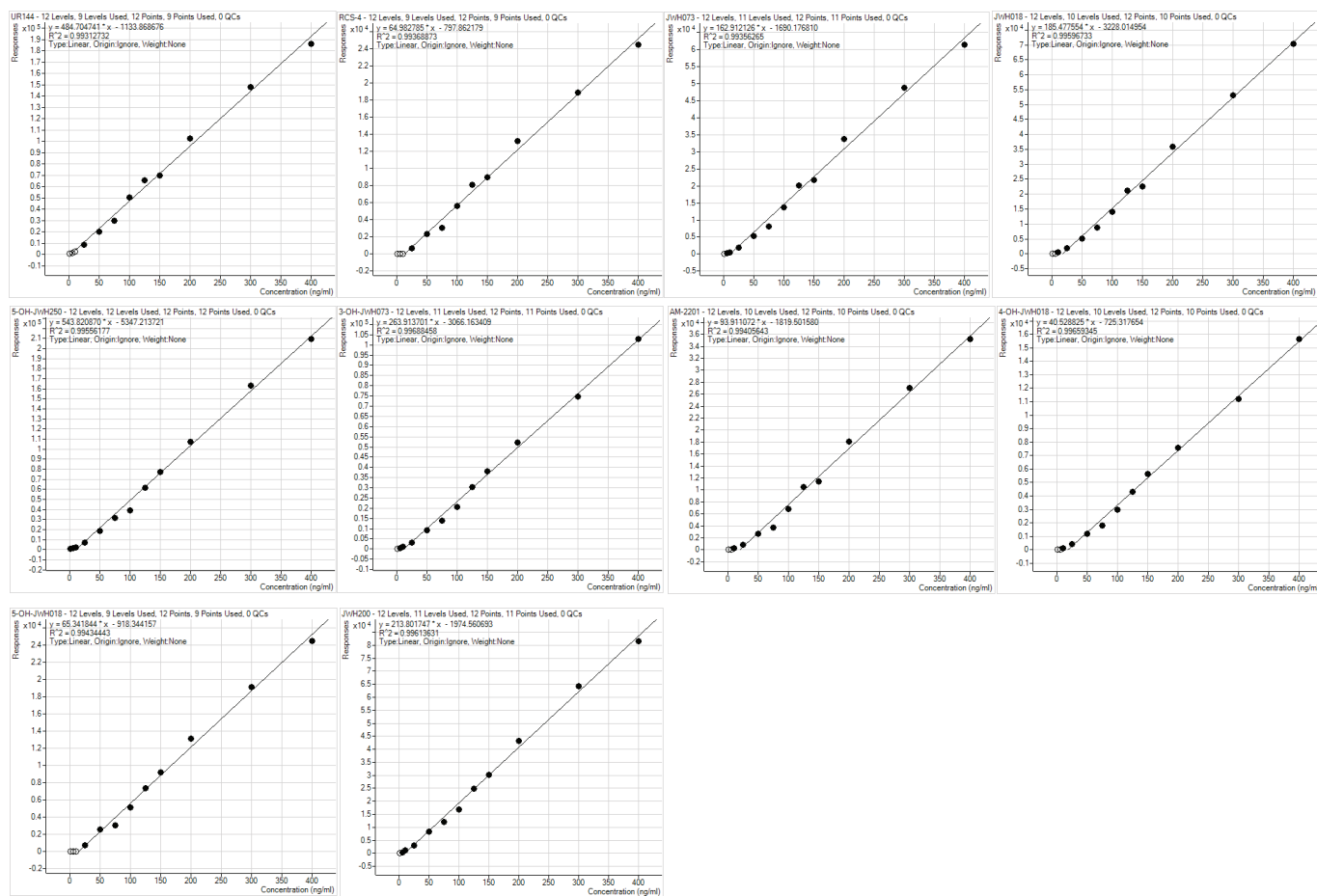

**Figure 2.** Calibration curves for extracted levels of spiked whole blood from 1 ng/mL to 400 ng/mL using 1 mL capacity ISOLUTE SLE+ columns showing R<sup>2</sup> values ranging from 0.9931 to 0.9968.

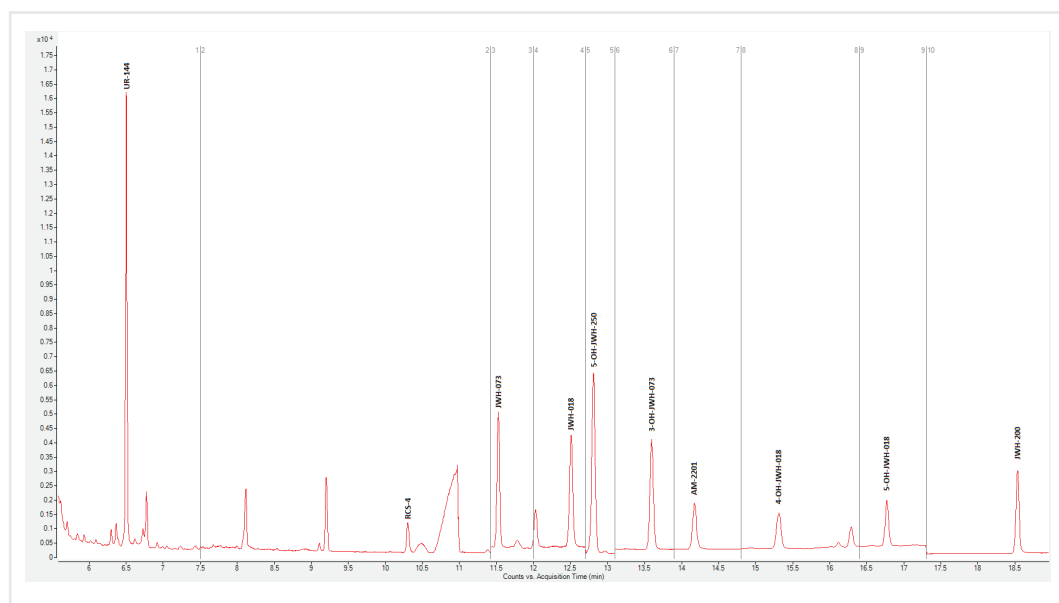

**Figure 4.** SIM chromatogram for whole blood spiked at 50 ng/mL

**Table 3.** Lower Limits of Quantitation (LLOQ) using the ISOLUTE SLE+ procedure

Analyte	Lower Limit Of Quantitation
UR-144	10 ng/mL
RCS-4	25 ng/mL
JWH-073	10 ng/mL
JWH-018	10 ng/mL
5-hydroxypentyl-JWH-250	25 ng/mL
3-hydroxybutyl-JWH-073	5 ng/mL
AM-2201	10 ng/mL
4-hydroxypentyl-JWH-018	10 ng/mL
5-hydroxypentyl-JWH-018	25 ng/mL
JWH-200	5 ng/mL

## Ordering Information

Part Number	Description	Quantity
820-0140-C	ISOLUTE® SLE+ 1 mL Sample Volume Columns	30
PPM-48	Biotage® PRESSURE+ 48 Positive Pressure Manifold 48 Position	1
SD-9600-DHS-EU	Biotage® SPE Dry Sample Concentrator System 220/240 V	1
SD-9600-DHS-NA	Biotage® SPE Dry Sample Concentrator System 100/120 V	1
C103198	TurboVap® 96 without racks 100/120 VAC	1
C103199	TurboVap® LV without racks 220/240 VAC	1

For the latest application notes visit [www.biotage.com](http://www.biotage.com)

### EUROPE

Main Office: +46 18 565900  
Toll Free: +800 18 565710  
Fax: +46 18 591922  
Order Tel: +46 18 565710  
Order Fax: +46 18 565705  
order@biotage.com  
Support Tel: +46 18 56 59 11  
Support Fax: +46 18 56 57 11  
eu-1-pointsupport@biotage.com

### NORTH & LATIN AMERICA

Main Office: +1 704 654 4900  
Toll Free: +1 800 446 4752  
Fax: +1 704 654 4917  
Order Tel: +1 704 654 4900  
Order Fax: +1 434 296 8217  
ordermailbox@biotage.com  
Support Tel: +1 800 446 4752  
Outside US: +1 704 654 4900  
us-1-pointsupport@biotage.com

### JAPAN

Tel: +81 3 5627 3123  
Fax: +81 3 5627 3121  
jp\_order@biotage.com  
jp-1-pointsupport@biotage.com

### CHINA

Tel: +86 21 2898 6655  
Fax: +86 21 2898 6153  
cn\_order@biotage.com  
cn-1-pointsupport@biotage.com

To locate a distributor,  
please visit our website at  
[www.biotage.com](http://www.biotage.com)

#### Part Number: AN773

© 2013 Biotage. All rights reserved. No material may be reproduced or published without the written permission of Biotage. Information in this document is subject to change without notice and does not represent any commitment from Biotage. E&OE. Product and company names mentioned herein may be trademarks or registered trademarks and/or service marks of their respective owners, and are used only for explanation and to the owners' benefit, without intent to infringe.

For more information visit [www.biotage.com](http://www.biotage.com).